

## ADDING A SECOND UNIT FOR Mettler Toledo JL Models

[www.balances.com](http://www.balances.com) is your source for Mettler Toledo Balances

The Mettler Toledo JL models can weigh in various modes. However, you can have only two units showing at a time. The scales come factory preset to weigh in grams only. To add a second weighing unit, follow the steps below.

Turn Scale On by pressing  $\rightarrow \frac{\text{O/T}}{\text{On/Off}} \leftarrow$  button. Display will read **0.0 g (or 0.00g)**

Press and HOLD **Menu** button until you see **MENU**. Release **Menu** button. Display will read **rESEt**.

Press a **Menu** button ONCE and release. Display will read **F none**.

Press a **Menu** again and release. Display will read **Std**

Press a **Menu** again and release. Display will read **Unit 1 g**

Press a **Menu** again and release. Display will read **Unit 2 g**

Now press **F button** (button with 2 arrows on it). The display will show **Unit 2 kg**

Again press **F button**. The display will show **Unit 2 ct**.

Repeat the last step until the display shows **Unit 2 and the desired weighing unit**.

STOP

Now Press and HOLD **Menu** button until you see **StorEd**. Release **Menu** button.  
The second unit has been added.

To switch between the two units, press **F button** (button with 2 arrows on it)

[www.balances.com](http://www.balances.com) is your source for Mettler Toledo Balances