

ADDING A SECOND UNIT FOR Mettler Toledo JL Models

www.balances.com is your source for Mettler Toledo Balances

The Mettler Toledo JL models can weigh in various modes. However, you can have only two units showing at a time. The scales come factory preset to weigh in grams only. To add a second weighing unit, follow the steps below.

Turn Scale On by pressing $\rightarrow \frac{\text{O/T}}{\text{On/Off}} \leftarrow$ button. Display will read **0.0 g (or 0.00g)**

Press and HOLD **Menu** button until you see **MENU**. Release **Menu** button. Display will read **rESEt**.

Press a **Menu** button ONCE and release. Display will read **F none**.

Press a **Menu** again and release. Display will read **Std**

Press a **Menu** again and release. Display will read **Unit 1 g**

Press a **Menu** again and release. Display will read **Unit 2 g**

Now press **F button** (button with 2 arrows on it). The display will show **Unit 2 kg**

Again press **F button**. The display will show **Unit 2 ct**.

Repeat the last step until the display shows **Unit 2 and the desired weighing unit**.

STOP

Now Press and HOLD **Menu** button until you see **StorEd**. Release **Menu** button.
The second unit has been added.

To switch between the two units, press **F button** (button with 2 arrows on it)

www.balances.com is your source for Mettler Toledo Balances